

# PRODUCT INFORMATION

## Formaldehyde Solidifier FSC-1

Fast absorption of formaldehyde spills

### Summary

- + **Easy to apply:** Open shaker, push paper protection, and generously sprinkle onto the spillage. Leave it to solidify for a short time, then absorb quickly with scoop and plastic scraper and close airtight in a waste bag / waste container with the cable tie. Thereafter, dispose according to the waste key 07 01 04\*.
- + **Application area:** Formaldehyde spills, 4% (e.g. pathology)
- + **Binding characteristics:** For Formaldehyde – No guarantee for other chemicals!
- + **Wear protective clothing when removing spills:**  
Wear tested PPE (gown, protective gloves, protective boots) of the highest category III with proved protection against Formaldehyde. Pay attention to the corresponding permeation times! In addition, you should wear a gas filtering respirator (half-mask or full face mask, no FFP masks) with appropriate filters and gas-tight goggles.
- + **Disposal:** Waste in need of supervision or need of special supervision (waste bag: 07 01 04\*); separately collect and dispose airtight!

### Version

Product	Formaldehyde solidifier FSC-1
Order number	100113

### Material characteristics

Absorbability	1 g FSC-1 = Ø 245 g Formaldehyde 4%
Shaker bottle – content	312 g solidifier

Shelf life	In closed container, dark and dry
------------	-----------------------------------

Producer/ Distributor	<b>Producer:</b> Safetec of America Inc., USA. <b>Distributor: Berner International GmbH,</b> Werner-von-Siemens-Str. 19, 25337 Elmshorn, Germany Tel: +49 4121 4356 - 0, Fax: +49 4121 4356 - 20, <a href="mailto:info@berner-safety.de">info@berner-safety.de</a> , <a href="http://www.berner-safety.de">www.berner-safety.de</a>
--------------------------	--