



Sterile, alcohol based ready to use solution for surface disinfection in cleanrooms.

perform® sterile alcohol IPA (WFI)

Our Plus

- fast, effective and residue free
- sterile (γ-irradiated) and microbially filtered (0.2 µm)
- suitable for cleanrooms grade A/B, ISO class 5/6
- easy to use in a handy, sterile 500 ml spray system
- with WFI (Water for Injection)
- colour and perfume free
- efficacy tested according to European norms (EN)

Application areas

Disinfection of surfaces and equipment / devices in cleanrooms, in particular for class A and B cleanroom areas, according to:

- Annex 1 of EU Guide to Good Manufacturing practice: Manufacture of sterile medicinal products, March 2009.
- FDA Guidance for Industry: Sterile drug products produced by aseptic processing - current Good Manufacturing practice; U.S. Department of Health and Human Services, Food and Drug Administration, September 2004.

Instructions for use

Apply perform® sterile alcohol IPA WFI undiluted to objects and surfaces and allow to take effect. Ensure complete wetting. The product dries off quickly. Maximum application rate 50 ml/sq.m. Do not use to treat surfaces that are particularly sensitive to alcohol, such as acrylic glass. perform® sterile alcohol IPA is also available with purified water (VE-water).

Microbiological efficacy

Efficacy	Concentration	Contact time
bactericidal EN1276, EN13697	ready-to-use	1 min.

Efficacy	Concentration	Contact time
yeastocidal EN1650, EN13697	ready-to-use	1 min.
virucidal against enveloped viruses <i>in accordance with DVV (German Registered Association for Combating Viral Diseases)//RKI Guideline</i>	ready-to-use	30 sec.



perform[®] sterile alcohol IPA (WFI)

Product data

Composition:

100 g solution contains the following active substances: 63.1 g Propan-2-ol (equates 70% (v/v)).

Special advice

Use biocides safely. Always read the label and product information before use.

For all of the related safety information please refer to the Safety Data Sheet. To request the Safety Data Sheet please E-mail: tech.uk@schuelke.com

The sterile 500 ml spray system offers the following advantages:

- convenient size for handling and storage
- ergonomically shaped spray pump for ease of use
- remains sterile after opening

The bottles are sterile (with irradiation indicator) and double wrapped individually. perform[®] sterile alcohol IPA comprises alcohol of pharmaceutical quality in accordance with the Ph. Eur.

Information for order

Item	Delivery form	Item no.
perform ster alc IPA WFI 500 ml FL	10/Carton	121101

Environmental information

schülke manufactures products economically and with advanced, safe and environmentally friendly production processes while at the same time maintaining out high quality standards.

Expert opinion and information

Please visit our website for an overview of all available literature/reports on all schulke products online at <http://www.schuelke.com/>

For any queries you might have about the product or using it please call 0114 2543500 or E-mail mail.uk@schuelke.com



Schülke & Mayr GmbH holds a Manufacturer's Authorisation according to sect 13 para 1 German Drug Law and Certificates of GMP Compliance for medicinal products.

schülke UK
Schülke & Mayr UK Ltd
Cygnet House, Meadowhall
Sheffield, S9 1AT
Phone +44-114-254 35 00
Fax +44-114-254 35 01
mail.uk@schuelke.com

schülke Headquarters
Schülke & Mayr GmbH
Robert-Koch-Str. 2
22851 Norderstedt, Germany
Phone +49 40 - 52100 - 0
Fax +49 40 - 52100 - 318
www.schuelke.com
info@schuelke.com